



## Coverslip Acid Wash

Fisher Microscope Cover Glass Circle

### Day 1

1. Place 30mm round glass cover slips in holders and immerse in 20%  $H_2SO_4$  overnight.

### Day 2

2. Remove from acid bath and immerse cover slips in holders in Milli-Q  $H_2O$  bath 5 minutes. Repeat this two times with fresh  $H_2O$ .
3. Remove the cover slips from  $H_2O$  bath, blot dry extra  $H_2O$  and place in oven at  $37^\circ C$  to dry.